

pH 10.00

Calibration solution

Works with any pH Probe

20ml each pouch

25 pouches per box

±0.02 pH at 25°C

**Standardized against
NIST-certified references**

2 year shelf life at room temperature



Pre calibration

Remove the soaker bottle and rinse off the pH probe. Once the probe has been thoroughly rinsed, dry off the probe with a paper towel.



Remove the top section of the pouch and insert your pH probe. Let the probe sit in the calibration solution until readings stabilize.

This process should take about 1 – 2 minutes.



1 – 2 mins

Once readings have stabilized you are now ready to calibrate.

Post calibration

Once the calibration is complete, dispose of the used solution.

